

# Magnetic Base

Date: 06/2020, v1.1

Adhesive back magnetic bases adhere to standard platens, allowing magnetic system cloths, abrasive discs or ferromagnetic discs to be utilized.

Size	Item Number
8" (203 mm)	90-208300
10" (254 mm)	90-208302
12" (305 mm)	90-208305



## Instructions

1. Clean the platen surface with acetone or isopropyl alcohol to remove any dirt or debris.
2. Remove the lining from the adhesive back and carefully apply the magnetic base to the plate. Press firmly to ensure that no bubbles form between the magnetic base and the platen surface.
3. The platen and magnetic base are now ready to use with any magnetic system cloths, abrasive discs or ferromagnetic discs.
4. To remove the magnetic base, lift the edge of the base until it completely separates from the platen.

## Maintenance

Periodically wipe the magnetic base clean to remove abrasive particles and suspensions. It can also be rinsed with water until it is clean; dry the base with clean air or a lint-free wipe.

Replace the magnetic base if it shows signs of wear, such as deep gouges, demagnetization, corrosion, nonuniformity due to overuse in one area or loss of adhesion between the base and platen.