

Glass Support Plate

Date: 06/2020, v1.2

The glass support plate is a smooth, clear 8" $\emptyset \times 0.25$ " T plate that provides an ultra-flat surface for lapping products (Figure 1). It is intended to be used on top of a standard or magnetic MultiPrepTM platen when standard platens do not provide a flat or smooth enough surface due to grooves and/or rough surfaces from manufacturing. Following standard alignment of the MultiPrepTM, the glass plate can be used accordingly:



Figure 1: Glass support plate (top), platen (bottom, not included with purchase)

Plain Glass Support Plate

- Place the glass plate on top of a dry standard MultiPrep™ platen.
- 2. Adhere the glass directly to the platen by wrapping electrical tape around the perimeter of both the glass and the platen (Figure 2).
- 3. Place diamond lapping film on the glass plate and use it as normal.



<u>Figure 2</u>: Electrical tape is wrapped around both the glass and the platen.

Magnetic Glass Support Plate

- Place a magnetic glass support plate on a magnetic MultiPrep™ platen (Figure 3).
- 2. Place diamond lapping film on the glass plate and use it as normal.
- 3. Tape and other adhesives are not necessary.

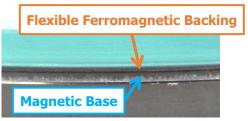


Figure 3: Magnetic glass plate on a magnetic platen

Lapping with a Glass Support Plate

The glass plate can be used as a lapping surface in conjunction with diamond compounds or diamond lapping films and a designated polishing lubricant. The plate is hard and does not compress, allowing for tighter edge tolerances and better flatness than compressible polishing cloths.

Product Numbers

69-10030 Glass Support Plate, Plain, 8"

69-10030M Glass Support Plate, Magnetic Back, 8"